

96-Well Vacuum Manifold

Catalog Number: 96Vac (96-Well Vacuum Manifold) \$375/each

Product Information:

The 96Vac, 96-Well Vacuum Manifold, is intended to be used with the vacuum protocols in our 96-Well Nucleic Acid Purification Kits. It can accommodate both the RNA and DNA 96-Well Plates.

The Lower Plate has a drilled hole on one side to accommodate the Vacuum Nozzle. Ensure that the Vacuum Nozzle has a Vacuum Nozzle Ring attached to its base, and screw it in until the Vacuum Nozzle Ring is firmly against the side of the Lower Plate. The Collection Chamber will fit in the open area in the center of the Lower Plate. It is easily removable, so that the user can discard flow-through during a procedure.

The Upper Plate fits securely on top of the Lower Plate, and each corner should be aligned, with the only separation between the two plates being made by the Lower Plate's sealing ring, located on its top.

The Upper Plate has a wider sealing ring, which will create a seal on the 96-Well Plates provided in a 96-Well Nucleic Acid Purification Kit.

Contents:

96Vac (96-Well Vacuum Manifold)	Quantity
Upper plate	1
Lower Plate	1
Collection Chamber	1
Vacuum Nozzle	1
Instruction Manual	1

Pictures for assembling the unit:

